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
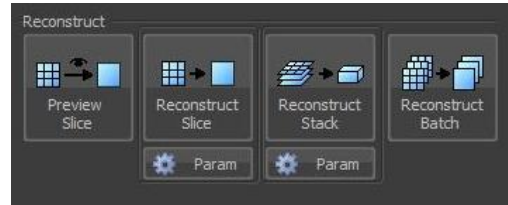
Number of steps and pitch sizes setttable on the N-SIM Z-Stack window (magnification: 100x)

Step Pitch	7	8 to 15	16 to 31	32 to 63	64 to 127	128 to 255	256 to 511
0.03 μm	No	No	Yes	Yes	Yes	Yes	Yes
0.06 μm	No	Yes	Yes	Yes	Yes	Yes	Yes
0.12 μm	Yes	Yes	Yes	Yes	Yes	Yes	Yes
0.24 μm	Yes	Yes	Yes	Yes	Yes	Yes	No
0.48 μm	Yes	Yes	Yes	Yes	Yes	No	No
0.96 μm	Yes	Yes	Yes	Yes	No	No	No
1.92 μm	Yes	Yes	Yes	No	No	No	No
3.84 μm	Yes	Yes	No	No	No	No	No
7.68 μm	Yes	No	No	No	No	No	No

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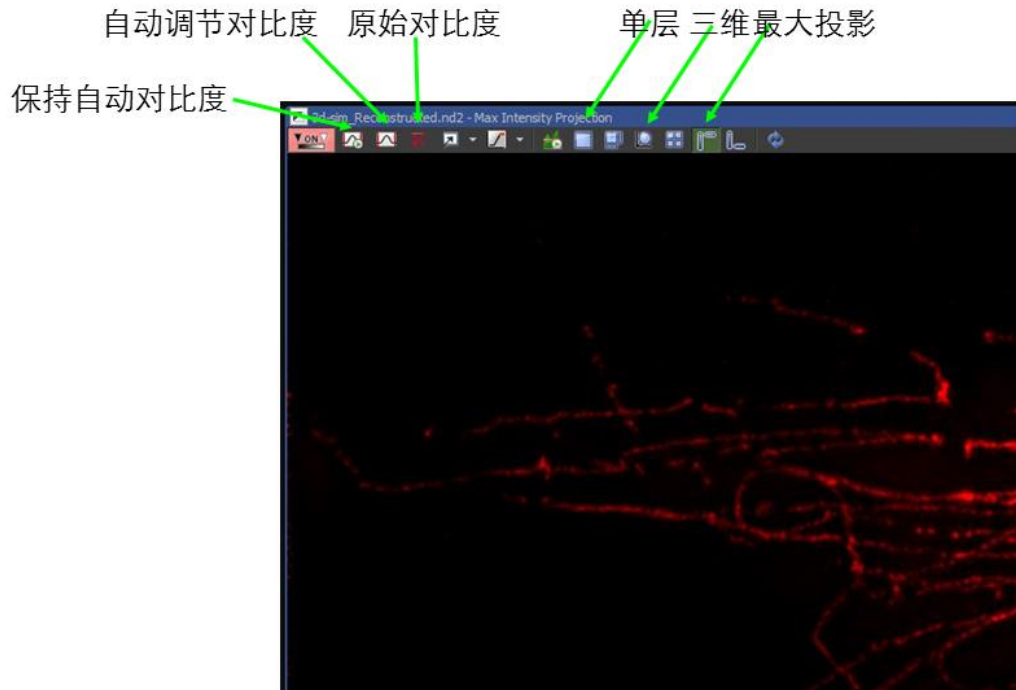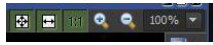
Methods for N-SIM:

3D-SIM images were acquired/recorded/performed on an N-SIM imaging system (Nikon) equipped with a 100X/1.49 NA oil-immersion objective (Nikon) and four laser beams (405, 488, 561 and 640). Images stacks with 0.12/0.24/0.48 μm interval were acquired and computationally reconstructed to generate super-resolution optical serial sections with two-fold extended resolution in both xy and z directions/in all axes. The reconstructed images were further processed for maximum-intensity projections and/or 3D-rendering with NIS-Elements AR 4.20.00 (Nikon).

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3D-SIM images were recorded using electron-multiplying CCD (charge-coupled device) cameras (Photometrics) on a DeltaVision OMX V3 imaging system (GE Healthcare) equipped with a $\times 100/1.40$ NA oil-immersion objective (Olympus) and three laser beams (405, 488 and 593 nm). Immunofluorescence images of Alexa Fluor 647 were acquired with deconvolution microscopy. To obtain optimal images, immersion oils with refractive indices of 1.512 and 1.520 were used for U2OS cells on glass coverslips and MTECs on Transwell polyester membrane inserts, respectively. The microscope was routinely calibrated with 100 nm fluorescent spheres to calculate both the lateral and axial limits of image resolution. Serial z-stack sectioning was carried out at 125 nm intervals. Images were reconstructed with the softWoRx 5.0 software package (GE Healthcare). The following parameters were set: channel-specific optical transfer functions, Wiener filter 0.001000, discard negative intensity background (post-processing), drift correction with respect to the first channel and custom K0 guess angles for camera positions. The pixel registration was corrected to be less than 1 pixel for all channels using 100 nm TetraSpeck beads (Life Technologies). The reconstructed images were further processed for maximum-intensity projections with softWoRx 5.0. 3D rendering was produced by using Imaris 7.6.0 (Bitplane).

All the antibodies used are listed in Supplementary Table 2.

Wide-field and structured illumination microscopy

Cells were fixed in methanol for 5 min at -20°C . Antibody incubations and washings were performed as described previously (Meraldi et al., 1999). Super-resolution 3D-SIM imaging was performed on a DeltaVision OMX V3 system (Applied Precision) equipped with a $100\times/1.40$ NA PlanApo oil immersion objective (Olympus), and 405, 488 and 592 nm diode lasers and Cascade II:512 EMCCD cameras (Photometrics). Exciting light was directed through a movable optical grating to generate a fine-striped interference pattern on the sample plane. Emitted fluorescence light was directed through a set of appropriate dichroic beam splitter and emission filter according to its wavelength range onto three cameras. Image stacks of typically $4\ \mu\text{m}$ height with 15 images per plane (5 phases, 3 angles) and a z-distance of $0.125\ \mu\text{m}$ were acquired and computationally reconstructed to generate super-resolution optical serial sections with two-fold extended resolution in all three axes. Color channels were carefully aligned using alignment parameter from control measurements with $0.5\ \mu\text{m}$ diameter multi-spectral fluorescent beads (Life technologies). Images were acquired in the sequential imaging mode to rule out bleed-through (supplementary material Fig. S1C). SI reconstruction and image processing was performed with the SoftWoRx 3.7 imaging software package (Applied Precision).

3D-SIM. NRK cells were plated on glass coverslips for 24 h and starved for 4 h. After immunostaining, 3D-images were acquired on an N-SIM using a $\times 100$ oil-immersion lens (numerical aperture 1.49) by Z-stack scanning at 0.2 or $0.4\ \mu\text{m}$ intervals. Laser lines at 488 and 561 nm were used for excitation using Nikon's N-SIM microscopy system, which can achieve a resolution of 100 nm along the x - y axis and 300 nm along the z axis. The instrument resolution is verified by 100 nm fluorescent spheres according to the Nikon procedure after installation.

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