

Axio Scan Z1 ZEN Blue Image Export

Axio Scan Z1 图像导出



Version 2013.11.13
This version is based on Zen blue 2012

Drafted by
ZHAO Yibing, AP Shanghai


目录



1 单张图像的导出

2 图像的批量导出


单张图像的导出

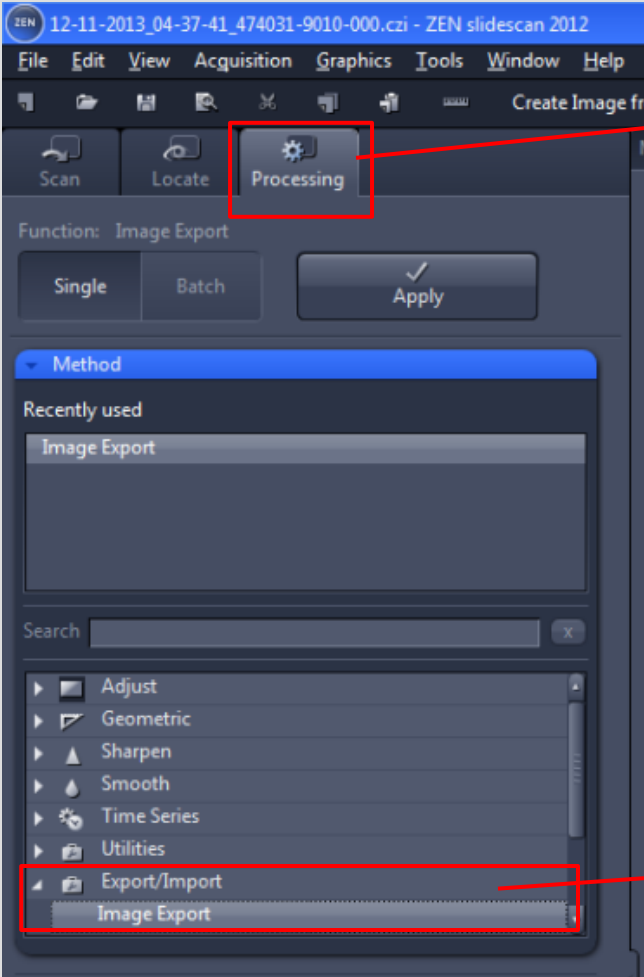


1. 选择选卡 “Processing”

2. 选择 “Export/Import”
→ “Image Export”


单张图像的导出

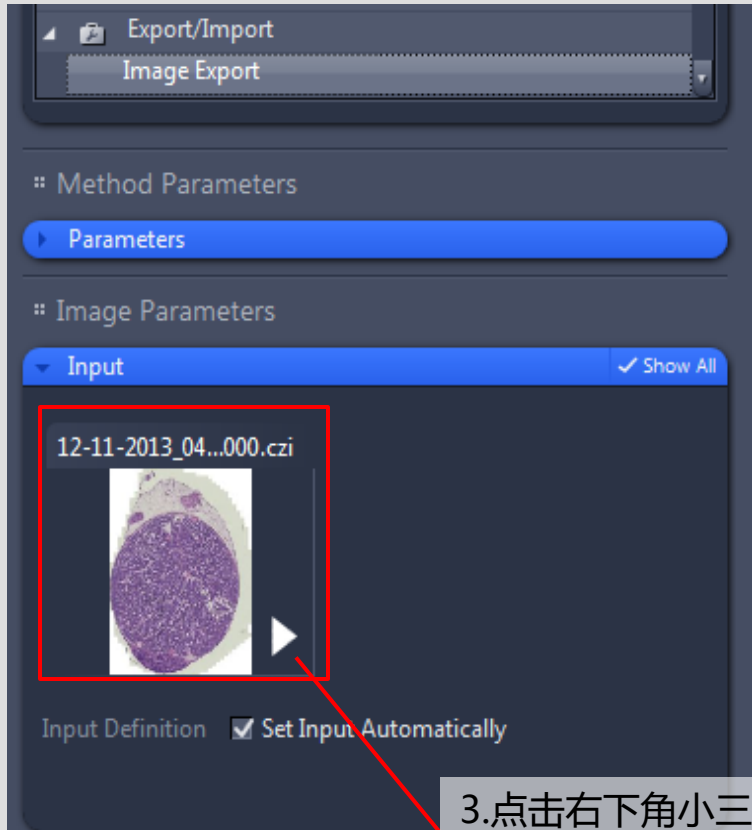

3. 点击右下角小三角，并选择需要导出的图像

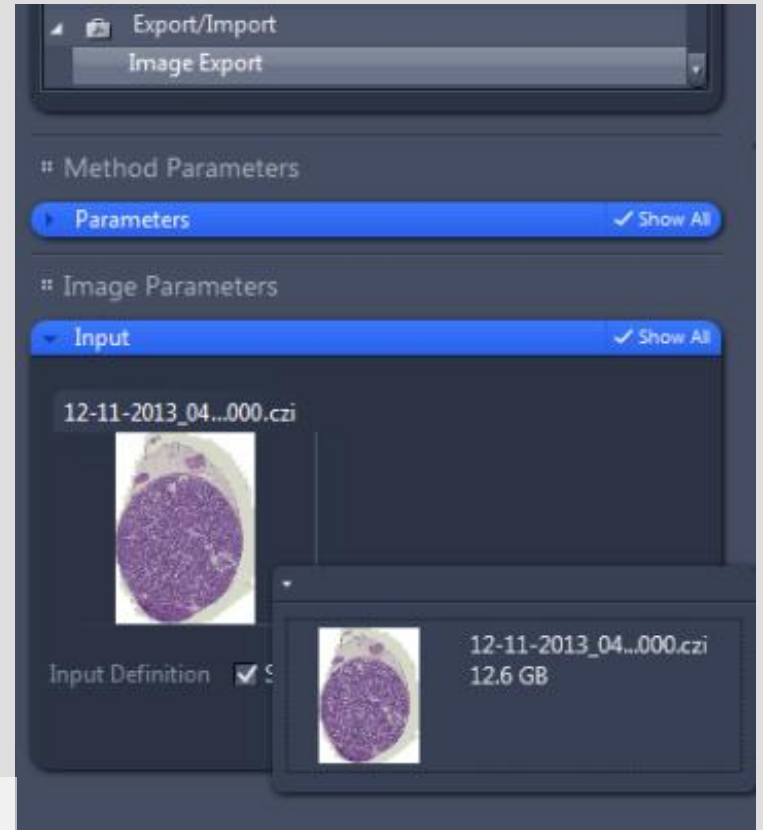

单张图像的导出

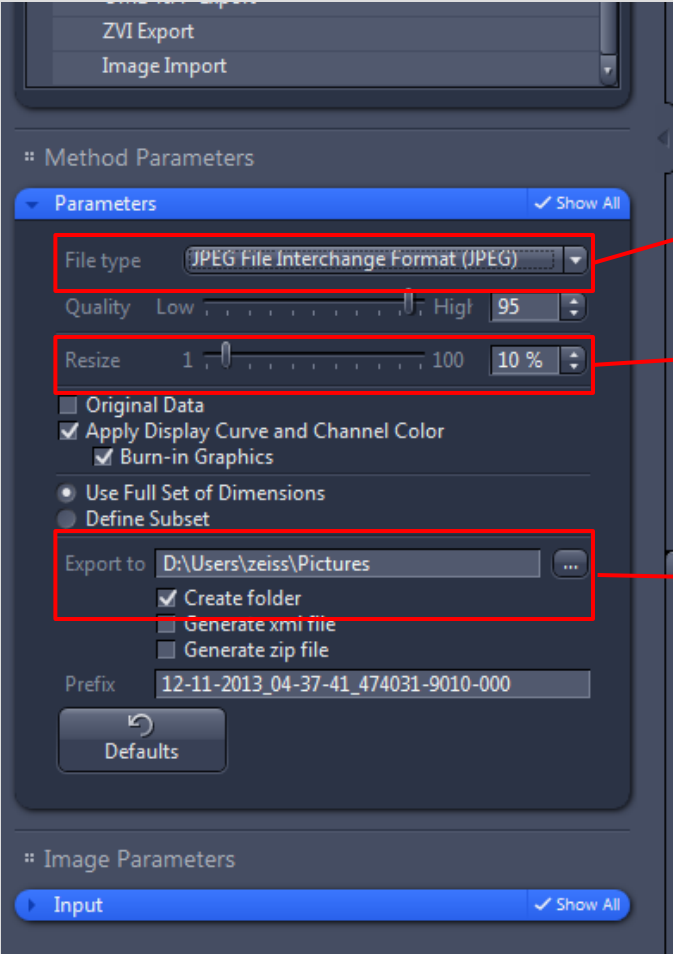

4. 选择需要导出的图像格式

5. 导出的图像大小，一般10%~20%

6. 选择保存的路径，可以选择是否需要建立新的文件夹

单张图像的导出



12-11-2013_04-37-41_474031-9010-000.czi - ZEN slidescan 2012

File Edit View Acquisition Graphics Tools Window Help

Workspace Zoom [Reset]

Design Dark Workspace Zeiss

Functions: Image Export

Single Batch **Apply**

Method

Recently used

Image Export

Search

Time Series

Utilities

Export/Import

Image Export

Movie Export

OME TIF-Export

ZVI Export

Image Import

Method Parameters

Parameters Show All

File type: JPEG File Interchange Format (JPEG)

Quality: Low High 95

Resize: 1 100 10 %

Original Data

Apply Display Curve and Channel Color

Summ Graphics

Use Full Set of Dimensions

Define Subset

Export to: D:\Users\zeiss\Pictures

Create folder

Generate xml file

Generate zip file

Prefix: 12-11-2013_04-37-41_474031-9010-000

Defaults

Image Parameters

Input Show All

2,000 µm

7. Apply !!

Dimensions Graphics

Zoom: 100% Auto Fit

Tools: Navigator Interpolation

Channels: Bright

Single Channel Range Indicator Quick Color Setup

Reuse

Display Show All

Histogram Spline Mode

Auto (Min/Max) Best Fit 2.00 0.01 Current Reset

Black: 0 Gamma: 1.00 0.45 1.0 White: 255

Scaling: 0.22 µm/pix (Calibrated) System Information: Free RAM 56.35 GB CPU 43 % Frame Rate: Pixel Value: Position: X: 32581 Y: 54466 Storage Folder: D:\Users\zeiss\Pictures User: zeiss

3:43 PM

图像的批量导出

1. 选择 "Batch"

2. 选择 "Image Export"

3. 选择需要导出的图像格式

4. 导出的图像大小，一般10%~20%

5. 选择保存在当前文件夹

6. 添加或移除需要导出的图像

7. Apply

S	Consistent	File Name	Size	Method	Output Name	Output Storage Path
?		G:\Training\2013.11.05-08 Si...	644.48 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\Carter
?		G:\Training\2013.11.05-08 Si...	37.79 MB	Change Scaling		G:\Training\2013.11.05-08 Singapore\Sample\Carter
?		G:\Training\2013.11.05-08 Si...	99.88 MB	Change Scaling		G:\Training\2013.11.05-08 Singapore\Sample\Slides Frank vor Auslieferung
?		G:\Training\2013.11.05-08 Si...	226.29 MB	Change Scaling		G:\Training\2013.11.05-08 Singapore\Sample\Slides Frank vor Auslieferung
?		G:\Training\2013.11.05-08 Si...	129.75 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder\Everytile
?		G:\Training\2013.11.05-08 Si...	131.01 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder\Number of Point
?		G:\Training\2013.11.05-08 Si...	1.37 GB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	1.08 GB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	1.34 GB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	129.61 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	15.23 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	1.61 GB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	1.7 GB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	12.53 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	104.49 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	176.21 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	323.17 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	340.42 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder
?		G:\Training\2013.11.05-08 Si...	148.33 MB	Image Export		G:\Training\2013.11.05-08 Singapore\Sample\New folder



We make it visible.